



Microfluidics™

analytik

Superior Knowledge | Superior Results



M110P

Microfluidizer®

Pharma Basic Processor

M110P Microfluidizer® Pharma Basic Processor

Microfluidizer Benchtop Processor Model is Recommended for:

- ◆ Production of stable nanodispersions and nanoemulsions
- ◆ Cell disruption (yeast, E. coli, etc.)
- ◆ Microencapsulation in polymers, liposomes & oils
- ◆ Deagglomeration

The M110P models have been designed to reliably achieve continuous operating pressures up to 30,000 psi. Microfluidizer™ processors maximize the energy-per-unit fluid volume, resulting in uniform submicron particles.



Microfluidics™

Standard Features

Interaction Chamber™ Material	Diamond
Plunger Material	Zirconia Ceramic
Enclosure	Stainless Steel
Drive Method	Electric/Hydraulic
Auxiliary Processing Module™	Ceramic
300 ml Glass Feed Reservoir	Yes
Cooling Coil Assembly	Yes

M110P “Plug & Play” Laboratory Model for Continuous High-Shear Processing

- ◆ Achieve processing pressures up to 2068 bar (30,000 psi)
- ◆ Produces product flow rate up to 120 ml/min (100 ml/min on 50Hz model)
- ◆ ASME BPE-compliant
- ◆ cGMP-ready
- ◆ Tube-in-tube pharma-grade heat exchanger
- ◆ 7” HMI color touchscreen control
- ◆ Inlet & outlet thermocouples with display on touchscreen
- ◆ Pressure transducer with display on touchscreen
- ◆ Highest shear rates on the market
- ◆ Guaranteed scalability
- ◆ Customizable options to suit every application
- ◆ CE-compliant

Utilizing Microfluidics’ fixed-geometry Interaction Chamber™ technology, and a ceramic (zirconia) plunger, the M110P Microfluidizer processor is capable of processing a wide variety of applications such as emulsions, suspensions, and cell disruption, including the most difficult yeast and plant cells, in as few as 1 - 2 passes. What’s more, the process is repeatable and is guaranteed to scale up to pilot and / or production volumes.

M110P Microfluidizer® Pharma Basic Processor

The M110P family of machines contains an on-board 1.5 kW (2 HP) electric-hydraulic drive that powers a single acting intensifier pump. Process pressures may be adjusted up to 30,000 psi (2068 bar) on the M110P to produce the desired product results.

As the reciprocating intensifier pump travels through its pressure stroke, it drives the product at constant pressure through the Interaction Chamber™. Within the Interaction Chambers™ are unique fixed-geometry microchannels through which the product stream will accelerate to high velocities. This creates high-shear and impact forces to bring about the desired results as the high velocity product stream impinges on itself and / or on wear-resistant surfaces.



Upon exiting the Interaction Chamber™, the product flows through an external heat exchanger, which precisely regulates the product to a desired temperature. At this point, the sanitary product may be recirculated through the system for further processing or sent to the next step in the process.

Specifications

Description	M110P Pharma Basic
Pressure Range	up to 30,000 psi (2068 bar)
Minimum Sample Size	50 ml
Flow Rate* Approximately (With Water)	Up To 120 ml/min – 60 Hz Up To 100 ml/min – 50 Hz
Product Temperature Limit	158°F (70°C)
Power Requirements	60 Hz/120 VAC/20 Amp 60 Hz/230 VAC/10 Amp 50 Hz/220 VAC/12 Amp
Dimensions L x W x H	40" x 27" x 23" (102 cm x 69 cm x 58 cm)
Weight	300 lbs (136kg)

Options

- Larger stainless steel reservoirs
- Seal quench system
- Additional Interaction Chamber™
- Additional Auxiliary Processing Module™ (APM™)
- Batch processing
- IQ/OQ Documentation and Execution
- FAT, SAT, on-site start-up & operator training
- Preventive maintenance service

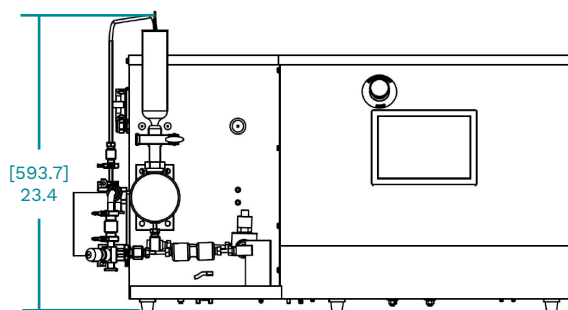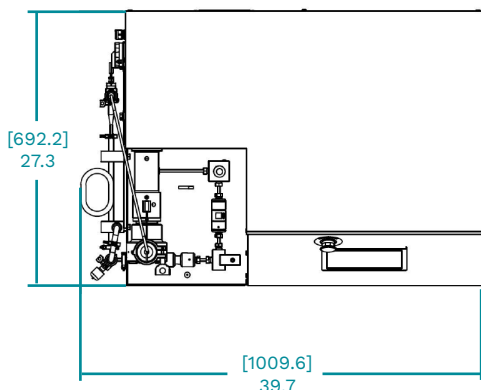

Manufactured by Microfluidics, distributed in the UK and Ireland by **analytik**.